



Electronic Crimpers and Decappers

Hand-held Rechargeable Crimpers and Decappers

- Precision, rechargeable crimpers and decappers for 8, 11, 13 & 20mm caps
- Easy to use, comfortable grip
- Hundreds of reproducible crimps from one charge
- Adjustable crimping force
- Includes standard Black & Decker VersaPak® Gold battery and charger

Reliable, Reproducible Crimping with the Push of a Button

Tired of crimping hundreds of vials with manual crimpers? Tired of losing your samples due to poorly crimped seals? The Electronic Crimper solves these problems and crimps autosampler crimp top vials with the touch of a button. Using the Black & Decker VersaPak® Gold rechargeable battery, this motor-driven, hand-held, high precision tool lets you crimp hundreds of vials per battery charge.

Designed to fit your hand, the electronic crimper and decapper allows a comfortable and natural position eliminating the fatiguing wrist and hand motion experienced when using manual, plier-type crimpers. The crimper jaws can be set electronically to the precise level required for reliable and reproducible crimping. The jaws are designed to fit around closely-spaced vials like those in standard vial holders or autosampler trays. The crimping force is set using up-and-down adjustment buttons and is controlled electronically with an internal microcontroller – giving you the reliability you need for a gas-tight seal on your samples.

The electronic crimper comes complete with the crimper, battery, and battery charger. A spare and/or replacement VersaPak® battery is also available.



Easy-to-Use

The comfortable grip makes the Electronic Crimper and Decapper easy-to-use. Simply place the jaws over the crimp cap and vial and push the cycle button with your gripping hand until the motor has engaged to cause a full crimping cycle. Within seconds, the user-adjustable jaws provide a precision crimp – a result that is repeated vial after vial. Works best when used in a seated position or standing at a GC autosampler tray.



- Size:** 13cm (L) x 4cm (W) x 18cm (H)
- Weight:** w/o battery: 445g, CE Certified w/battery 595g, charger: 290g
- Body:** Glass-filled nylon
- Jaws:** Carbon steel, electroless nickel plated
- Battery:** Black & Decker VersaPak Gold NiMH
- Voltage:** 3.6 V, Full charge time: 6-9 hrs.

Description	Cat. No.
Electronic Crimper, 8mm, Complete Kit *	16-100000
Electronic Crimper, 11mm, Complete Kit *	16-100002
Electronic Crimper, 13mm, Complete Kit *	16-100004
Electronic Crimper, 20mm, Complete Kit *	16-100006
Electronic Decapper, 11mm, Complete Kit *	16-100008
Electronic Decapper, 13mm, Complete Kit *	16-100010
Electronic Decapper, 20mm, Complete Kit *	16-100012

Accessories

Replacement VersaPak® Gold (NMH) Battery	16-100014
--	-----------

* State country of use when ordering



Greyhound Chromatography & Allied Chemicals
 6 Kelvin Park, Birkenhead, Merseyside, CH41 1LT, UK
 Tel: (+44) 0 151 649 4000 • Fax: (+44) 0 151 649 4001
 Email: info@greyhoundchrom.com • Web: www.greyhoundchrom.com